

IMMUNOSTAINING BY HRP AND AP Protocols

STAINING BY PEROXIDASE (IN SOLUTION)

Detection of peroxidase activity (Immunoassay):

Dissolve 1 mg TMB (Art. No. 6350) in 0.1 ml dimethylsulfoxide (Art. No. 4720).
Add 9.9 ml of a 0.1 M sodium acetate solution (pH 6.0) (Powder: Art. No. 6779), filter
and add H_2O_2 (Art. No. 8070) (final concentration 0.01 %).

Always prepare freshly!

Incubate 10-30 mins at room temperature (approx. 50 μl per microtitre well)
finally add 50 μl 1 M H_2SO_4 , Art. No. X873, per well
Photometric quantitation at 450 nm.

Reference: Bos, E.S. *et al.*, (1981) *J. Immunoassay* 2:187.

STAINING BY ALKALIC PHOSPHATASE (PRECIPITATE)

Assay protocol for use of BCIP/NBT in immunoblot procedures:

Stock solutions (all solutions are stable for at least 1 year at 4 °C):
0.5 g NBT (Art. No. 4421) in 10 ml 70 % dimethylformamide (Art. No. T921);
0.5 g BCIP p-toluidine salt (Art. No. 6368) in 10 ml 100 % dimethylformamide.
Incubation buffer for alkaline phosphatase:
100 mM NaCl, 5 mM MgCl_2 , 100 mM Tris (pH 9.5).

Fresh substrate solution:

66 μl NBT stock solution + 10 ml incubation buffer, mix well,
add 33 μl BCIP stock solution. Use within 1 hour.

Blot development:

Approx. 10 ml substrate solution per 15 x 15 cm^2 membrane surface.
Develop at room temperature until bands become visible (approx. 30 mins).


Reaction stop:

Rinse with PBS/20 mM EDTA.


NBT / BCIP Reaction Mechanisms

Alkaline phosphatase mediated reaction of 5-bromo-4-chloro-3-indolylphosphate (BCIP or X-Phosphate) and Nitroblue tetrazolium salt (NBT). By enzymatic dephosphorylation, BCIP is converted to indoxyl which subsequently tautomerizes to the ketone. In basic pH values, this ketone dimerizes, resulting in a weak blue colour. During dimerization, the ketone also releases H^+ ions, which reduce the colour enhancer NBT to purple-coloured diformazan. Both dyes precipitate in the direct vicinity of the phosphatase, hence staining this site in a dark violet.


Reaction:

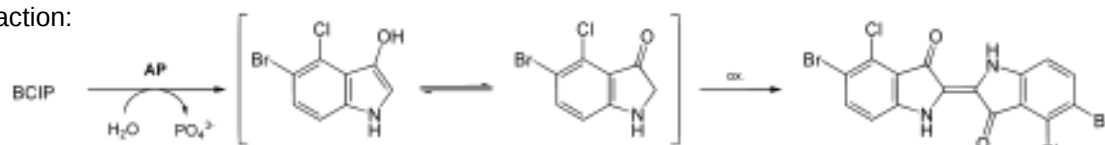

BCIP:

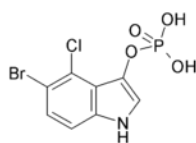

NBT:

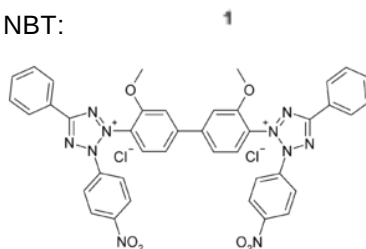

Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5
76185 Karlsruhe
Postfach 100121
76231 Karlsruhe
Telefon: +49 (0) 721/ 5606-0
Telefax: +49 (0) 721/ 5606-149
E-Mail: info@carlroth.com
Internet: www.carlroth.com

s.s. 03/2016